

## PIPETBOY *comfort*



The new superior **PIPETBOY** with charging station  
Outstandingly flexible, lightweight and durable!

# PIPETBOY *comfort*

## Light as a feather

A lightweight 190 grams, the strain is taken out of even prolonged pipetting work. Hands, arms and shoulders remain fatigue-free longer, enhancing your comfort.



## Charger and stand – a

Either with or without a pipette, utmost flexibility and ease of use. **PIPETBOY *comfort*** is on standby when you need it.


## Ergonomically designed

The **PIPETBOY *comfort*** sits snugly in the hand naturally matching its shape. The rounded handle also fits perfectly in small hands. Pipetting has never been so comfortable.



## Flexible and compatible

The silicone Pipette Mount guarantees a sure grip on practically all available plastic or glass pipettes, ideally in the 1 – 50 ml volume range.

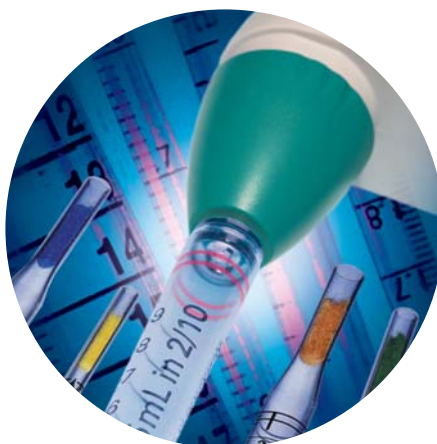

## Maximum Safety

The easy-to-change sterile filter provides your **PIPETBOY *comfort*** with reliable protection from liquid contamination and damage. And if you do happen to aspirate critical liquids with the pipette, the **PIPETBOY *comfort*** silicone Pipette Mount, the Pipettor Nose Piece and the filter rubber are all autoclavable while the sterile filter can be replaced. So easy and so safe.





classic



volcano




pacific



sahara


## All in one

on the table or on the wall – the  
of use for your laboratory. Your  
and fully charged, whenever you

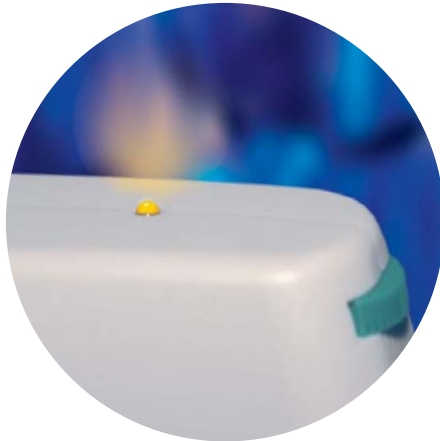

## Battery indicator

An illuminated yellow LED shows when the battery needs to be charged. While the battery is being charged, the LED flashes to show that the contact has been correctly established. You know the status of your rechargeable battery at all times. When the **PIPETBOY comfort** is connected directly to the mains adapter, you can still pipette while it is charging. The **PIPETBOY comfort** is always on standby!

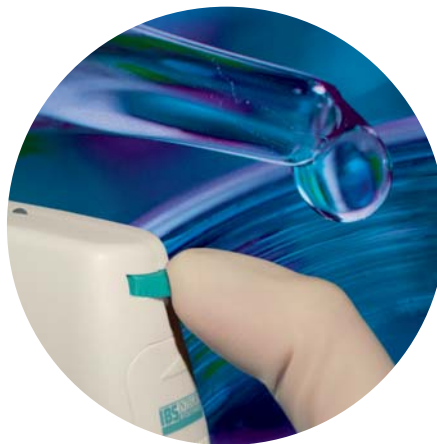

## From rapid pipetting to sensitive precision dosing

You can control the speed of pipetting continuously via the knurled screw. Precise dosing is also possible by applying different amounts of pressure to the triggers.



## Powerful

2 NiMH rechargeable batteries deliver optimum performance ensuring maximum power and durability with minimal weight. Fully charged batteries provide 8 hours of continuous operation – a truly masterly performance in view of the light weight. The **PIPETBOY comfort** can either be connected directly to the mains adapter or recharged via the charging console.



## Held perfectly in all positions

Whether on the wall or tabletop, inside or outside a hood, the **PIPETBOY comfort** is held perfectly – with or without a pipette.


**Wall mounting:** Magnet, adhesive strip or screw – the charger can be easily and securely mounted on any wall. The aerodynamic shape of the charger ensures minimal air turbulence in the laminar flow and reduces the risk of contamination. **Tabletop mounting:** Simply place the charger stand of your on top of your table – 4 rubber feet ensure stability.

## Ordering information

Description	Unit	Article No.
<b>PIPETBOY comfort classic</b> , incl. charging station, adapter, sterile filter, charger mounting kit, battery, one original Swiss knife	1	156 000
<b>PIPETBOY comfort volcano</b> , as Art. No. 156 000	1	156 001
<b>PIPETBOY comfort pacific</b> , as Art. No. 156 000	1	156 002
<b>PIPETBOY comfort sahara</b> , as Art. No. 156 000	1	156 003

## Ordering information for single and spare parts

No.	Description	Unit	Article No.
1	Filter rubber	1	153 225
2	Filter sterile, 0.45 µm	1	153 015
	Filter unsterile, 0.45 µm	1	153 016
3	Pipette Mount, silicone	1	151 020
4	Pipettor Nose Piece, (classic)	1	156 100
	Pipettor Nose Piece, (volcano)	1	156 101
	Pipettor Nose Piece, (pacific)	1	156 102
	Pipettor Nose Piece, (sahara)	1	156 103
5	Battery-Set, 2 Pieces (1.2 V AAA NiMH / 700 mAh)	2	156 151
6	Battery cover Set (classic)	1	156 110
	Battery cover Set (volcano)	1	156 111
	Battery cover Set (pacific)	1	156 112
	Battery cover Set (sahara)	1	156 113
7	Charging station (classic)	1	156 050
	Charging station (volcano)	1	156 051
	Charging station (pacific)	1	156 052
	Charging station (sahara)	1	156 053
8	Charger mounting kit (tape, magnet, base pads, screw)	1	156 150
9	Adapter:		
	USA, 120 VAC	1	156 630
	Europe, 230 VAC	1	156 631
	UK, 240 VAC	1	156 632
	Australia, 240 VAC	1	156 633
	Japan, 100 VAC	1	156 634

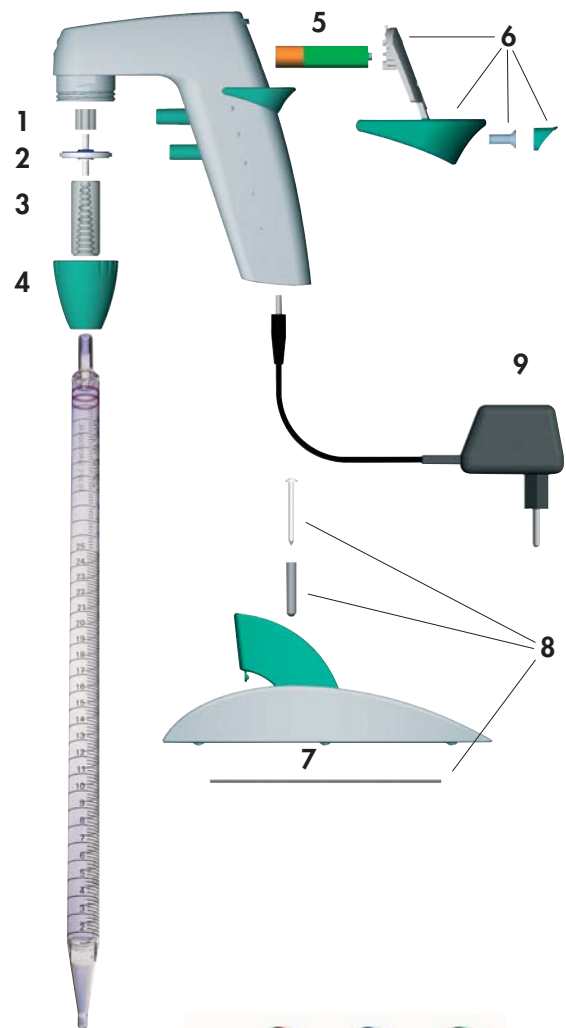

## Technical Data

Dimensions (LxHxW):	<b>PIPETBOY comfort:</b>	130 x 145 x 35 mm
	Charging station:	175 x 75 x 80 mm
	Mains adapter:	75 x 45 x 45 mm
Weight:	<b>PIPETBOY comfort:</b>	190 g
	Charging station:	240 g
	Mains adapter:	190 – 300 g
Batteries:	Rechargeable batteries	
		2 x 1.2 V AAA NiMH / 700 mAh
Operating condition:	Operation:	5 – 40° C, max. 80% RH
	Storage:	-10 – 50° C, max. 95% RH

AC adapter:	Input voltage:
USA 120 VAC	120 V~ / 60 Hz / 52mA / 6 VA
Europe 230 VAC	230 V~ / 50 Hz / 26mA / 6 VA
UK 240 VAC	240 V~ / 50 Hz / 26mA / 6 VA
Australia 240 VAC	240 V~ / 50 Hz / 26mA / 6 VA
Japan 100 VAC	100 V~ / 50 Hz / 60mA / 6 VA
	Output voltage:
all	9 V ~ / 300 mA / 2.7 VA

## INTEGRA Biosciences AG

CH-7000 Chur, Switzerland  
 Phone: +41 81 286 95 30  
 Fax: +41 81 286 95 33  
 E-mail: info@integra-biosciences.com  
 www.integra-biosciences.com

